

Name: Fusaric Acid

Alt. Name: Maloy's

Type:

<u>Components</u>	<u>Amount</u>
Agar	15g
Tryptone	5g
Yeast Extract	5g
Chlortetracycline *	.05g
NaCl	10g
NaH ₂ PO ₄ H ₂ O	10g

Method of Preparation:

*(12.5 mg / ml) Autoclave 20 min., cool to 45 °C, and then add 5 ml 20mM ZnCl₂ and 6 ml 2mg/ml fusaric acid. Can weigh out and autoclave with others or add as liquid from stock solution (stored in dark at 4 °)

Reference:

Source: