

Name: Green Indicator Plates

Alt. Name: Bresch's recipe modified 3-86

Type:

<u>Components</u>	<u>Amount</u>
Tryptone	6g
Nutrient Agar	20g
Yeast Extract	10g
NaCL	2g
Glucose	12g
H ₂ O	1000ml

Method of Preparation:

Autoclave; cool to 50 °C. Add the following components just before pouring plates: 4ml of 20%(wt/vol) sol'n Na₂CO₃* 10H₂O , 10ml of .5% water soluble blue, and 20ml of 2.5% Superchrome Yellow. Pour medium thickness.

Reference:

Source: